








# CONSOLIDATED STERILIZER SYSTEMS

## Medium Lab Series Steam Sterilizers General Specifications

Models		Temperature
<b>SR-24C</b> (24" x 36" x 36")		212°F-270°F
<b>SR-24D</b> (24" x 36" x 48")		212°F-270°F
<b>SR-24E</b> (24" x 36" x 60")		212°F-270°F
<b>SR-24F</b> (24" x 36" x 72")		212°F-270°F
<b>SR-26B</b> (26" x 26" x 49")		212°F-275°F

### General Specification

Steam Sterilizer, Radial-Arm Door(s), Hinged,  
Single Chamber, Double Wall

Choose from a variety of sizes and programmable control options for pre-vacuum or gravity operation. Consolidated sterilizers offer a range of performance options to meet the most demanding applications in clinical, animal and life science, biotechnology, pharmaceutical, and commercial/industrial applications.

*\* Exact upper temperature limit is model specific and may vary depending on model and/or options. Contact Consolidated for a solution tailored to your requirements.*

### Table of Contents

- Model Sizes and Weights..... 2
- Sterilizer Construction..... 2
- Sterilization Cycles..... 6
- Options..... 8
- Accessories..... 9
- Validation and Installation..... 10
- Preventative Maintenance..... 10
- Footprint Drawings..... 11
- Utility Information..... 13

### Features and Benefits

#### Simplified Maintenance, Low Cost of Ownership.

All Consolidated sterilizers are manufactured in the USA and built from commonly available parts to allow quick and cost effective field-level service and maintenance.

#### Serviceability.

Easy access to replaceable components, local component availability and common electrical and plumbing parts permit qualified facility or area service companies to maintain the sterilizer.

#### Control Flexibility.

A choice of programmable controllers allows a broad range of performance functions, complete with alarm, monitoring and communications required for internal or third-party compliance.

#### Performance Cycles—Basic to Advanced.

The fully-jacketed sterilizer design permits vacuum and pressure control when configured for pre-vacuum, post-vacuum, and more sophisticated functions such as air-over-pressure. Consolidated sterilizers are ideal for sterilizing wrapped and unwrapped goods, liquids, waste, and other applications.

#### Green and Environmentally Friendly.

Unique, new technologies reduce water and energy consumption without compromising performance.








Consolidated Medium Lab Series Sterilizers are available in single door, and pass-thru models. A versatile control system offers a range of performance options to meet the most demanding applications in life science, biotechnology, pharmaceutical, and commercial/industrial applications. Model SR-24D-X1 shown with X1™ control system.






## MODEL SIZES & WEIGHT

### Summary Selection Chart

Consolidated Medium Lab Series sterilizers are based on six chamber sizes, both single door and pass-thru models. Left or right door hinge side must be specified when ordering.

Model	Chamber Dimensions (w x h x f-b)	Volume (cu. ft. / liters)	Ship Weight (lbs/kg)	Ship Dimensions (w x h x f-b)
Single Chamber Models				
    				
<b>SR-24C</b>	24" x 36" x 36" 61 x 91.4 x 91.4 cm	18 cu. ft. 510 liters	2,700 lbs 1,228 kg	46" x 78" x 65" 116.8 x 198.1 x 165.1 cm
<b>SR-24D</b>	24" x 36" x 48" 61 x 91.4 x 122 cm	24 cu. ft. 680 liters	3,150 lbs 1,432 kg	46" x 78" x 78" 116.8 x 198.1 x 198.1 cm
<b>SR-24E</b>	24" x 36" x 60" 61 x 91.4 x 152.4 cm	30 cu. ft. 850 liters	3,600 lbs 1,637 kg	46" x 78" x 89" 116.8 x 198.1 x 226.1 cm
<b>SR-24F</b>	24" x 36" x 72" 61 x 91.4 x 183 cm	36 cu. ft. 1,019 liters	4,300 lbs 1,955 kg	46" x 78" x 98" 116.8 x 198.1 x 248.9 cm
<b>SR-26B</b>	26" x 26" x 49" 66 x 66 x 124.5 cm	19.2 cu. ft. 544 liters	2,700 lbs 1,227 kg	46" x 77" x 79" 116.8 x 195.6 x 201 cm

### Single Chamber, Pass-thru Models

    				
<b>PT-SR-24C</b>	24" x 36" x 36" 61 x 91.4 x 91.4 cm	18 cu. ft. 510 liters	3,150 lbs 1,432 kg	46" x 78" x 74" 116.8 x 198.1 x 188 cm
<b>PT-SR-24D</b>	24" x 36" x 48" 61 x 91.4 x 122 cm	24 cu. ft. 680 liters	3,600 lbs 1,637 kg	48" x 78" x 87" 122 x 198.1 x 221 cm
<b>PT-SR-24E</b>	24" x 36" x 60" 61 x 91.4 x 152.4 cm	30 cu. ft. 850 liters	4,050 lbs 1,841 kg	48" x 81" x 98" 121.9 x 205.7 x 248.9 cm
<b>PT-SR-24F</b>	24" x 36" x 72" 61 x 91.4 x 183 cm	36 cu. ft. 1,019 liters	4,750 lbs 2,160 kg	48" x 81" x 104" 122 x 205.7 x 264.2 cm
<b>PT-SR-26B</b>	26" x 26" x 49" 66 x 66 x 124.5 cm	19.2 cu. ft. 544 liters	3,050 lbs 1,386 kg	46" x 77" x 79" 116.8 x 195.6 x 201 cm

### High Performance Sterilization

Consolidated sterilizers are used in a wide range of applications including clinical, animal and life science, biotechnology, pharmaceutical, and commercial/industrial markets. Consolidated sterilizers are suitable for conventional or pre-vacuum sterilization for quality control, safety and regulatory compliance. In addition, all units can be customized to suit any unique requirements and applications of a particular facility.

Consolidated uses only commonly-available, open-source parts in all phases of sterilizer design and construction, thereby permitting quick, economical and locally-managed service throughout the life of the product. With maximum uptime engineered into the product, Consolidated sterilizers offer a superior return on investment as demonstrated through thousands of installations worldwide.

## STERILIZER CONSTRUCTION

### Sterilizer Chamber

Consolidated steam sterilizer vessels are manufactured with an inner chamber and a full outer jacket. The inner chamber is available in either type 316L stainless steel or nickel-clad steel construction. Jacket construction is available in carbon steel, type 304L stainless steel or type 316L stainless steel. Material selection is directly related to the chemical properties of available water supply and intended application. All vessels with type 316L stainless steel inner chambers are polished to a high luster finish and passivated. Sterilizers are equipped standard with a 1/2" NPT port for use with validation or process measurement. Pressure vessels are welded in compliance with ASME Section VIII, Division I.

### Chamber Door and Gasket System

Consolidated's sterilizer doors are designed with redundant, independent mechanical and software features to ensure maximum safety. Door design specifications meet all ASME code requirements.

- Consolidated hinged door autoclaves use solid silicone gaskets that do not require high pressure air, steam or vacuum to operate. Instead, the act of closing the door compresses the silicone gasket to create a secure and reliable seal. In the event that the gasket fails, the door can still be easily opened and the gasket quickly replaced.
- Multiple door arms assure that, in the event of a failure of any one arm, the door will continue to support the load within performance specifications.
- Door engagement is clearly visible to confirm closure.
- An electrical door switch ensures that the door is fully closed and locked prior to the start of the cycle. If the signal is lost during a sterilizing cycle, the cycle is automatically aborted.
- The door pressure clutch assembly is activated by chamber pressure to lock the door. Chamber pressure forces a membrane against the clutch plate, engaging the lock to permit tightening, but not loosening of the door.
- As the handle is turned to the closed position, the latching arms extend out into the door end-frame. Full rotation of the door handle presses the door against the chamber sealing surface thereby compressing the seal, and activating the electrical door limit switch.
- Pass-thru (double-door) sterilizers may be ordered with optional door interlocks. This electro-mechanical lock on each door minimizes the chance of cross-contamination between contained and uncontained areas by preventing both doors from being opened simultaneously. Additionally, this feature prevents the uncontained side door from opening until a sterilization cycle has successfully completed.

### Components

- Piping, fittings, and valves are available in stainless steel or brass/bronze. All parts and components subject to repair are publicly available (i.e. non-proprietary) from open market sources, including controls, valves and fittings.
- Safety relief valves and drains are integral to the plumbing configuration.
- Baffled steam inlets minimize direct condensate on the load, reduce load wetting, and help assure proper steam uniformity within the chamber.
- Cabinet insulation retains heat and minimizes heat loss (see Table 10) to the room. The sterilizer jacket and all steam service piping include a minimum 1" fiberglass-based insulation.

- On units equipped with steam generators, a manual reset pressure switch is included to back up the safety relief valve by automatically shutting down the steam generator if an overpressure condition exists.
- Easy-to-read pressure gauges ensure simple and safe operation.
- Other components required for the intended application may include a steam generator, vacuum pump, water ejector, etc. depending on the options chosen.

### Sterilizer Mounting

Consolidated sterilizers are available for installation as floor-standing cabinet-enclosed units with a stainless-steel superstructure to conceal plumbing and wiring. These sterilizers are also available for installation as recessed into one or two walls. Hook-ups for utility connections can be located anywhere on the unit as required, however house connections must be within 5 feet of the unit unless otherwise specified prior to installation.

For Biological Safety Level 3 installations a Bioseal® can be installed to maintain the contamination integrity of the clean side. Utilities can be mounted at the non-contained side of the installation and in the front, rear or designated side of the cabinet.

### X1™ Control and Monitoring System

Consolidated offers a state-of-the-art control system built from industry standard, open source components for international availability of parts and service. Consolidated's X1™ controllers offer proven reliability, based on an industrial PLC platform programmed specifically for Consolidated's sterilizers, complete with modern user interface and an intuitive touchscreen display (see chart on page5).

- Permits industry standard and custom sterilizer cycles.
- Simplifies operation with intuitive operator interfaces. Optional features allow the operator to add favorite shortcuts for frequently-used cycles and prevent users from modifying sterilization parameters on individual cycles.
- Assures quality control and process integrity.
- Features sterilizer cycle quality reporting and optional printed documentation.
- Internal battery retains all data in memory for at least 6 months.

### Alarms

Detailed performance and function alarms are included with the X1™ controller, as well an alarm history report visible on the touchscreen. Standard safety alarms are included in all Consolidated sterilizers. Process alarms and errors are documented and communicated both audibly and visually.

Service and Preventative Maintenance Reminder Alerts will be displayed on the touch screen and printed at adjustable use-based or time-based intervals. If desired, this functionality can be disabled.

### Cycle Safeguards

The following cycle safeguards help assure safe operation.

- No cycle can start unless the door is properly closed and locked.
- Control inputs automatically reject incorrect cycle parameters.
- Chamber gauges are installed on the operating end of the sterilizer to display pressure in psig and vacuum in inches-Hg.
- Exhaust is automatically condensed and cooled to a safe drainage temperature.
- Multi-rate liquid exhaust allows users to select an exhaust rate above and below an adjustable temperature set-point. This prevents boil-over of liquids and reduces total cycle time.
- Emergency Stop button cuts power to all 120VAC outputs and stops all processes.

### Calibration

All temperature and pressure sensors on the sterilizer can be calibrated using an intuitive three-point calibration method.

### Printer

A thermal printer is integrated into the main control panel. Printout includes all important information regarding the sterilization cycle recorded at user-defined intervals. Information includes: operator identification, cycle type, cycle parameters, cycle phase, time, chamber pressure and temperature, alarms or messages and cycle completion status.


### Electronic Data Storage via USB

Cycle data is stored on a non-volatile flash memory device and saved in CSV format. The flash memory device can be used to transfer the data to other computers or tablets. The user can use the data to create graphs and perform data analysis. This feature allows users to extract Raw Cycle Data (Time, Temp, Pressure), Cycle History (User ID, Cycle Name, Time, Temp, Alarm Code, Cycle Result), and Duplicate Cycle Data (an electronic copy of the paper print-out).

### Sustainable Features (Included with Every Autoclave)

#### Water Conservation & Monitoring System

All units are equipped with an integral WaterEco® water saving system that cools effluent discharge temperature (to drain) to less than 140° F (60° C) while keeping water consumption to a minimum (reduces cooling water by up to 90% over conventional sterilizers). This system also monitors, measures and reports quenching-water consumption through the touch screen.

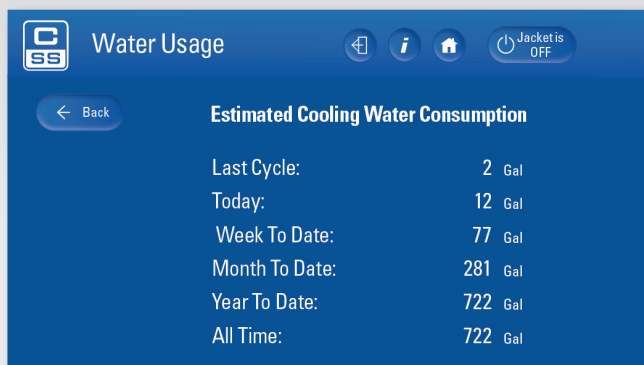

X1™ User-Interface: Water Usage Screen

#### EcoCalendar™

The X1 controls are equipped with a calendar-based automatic start-up and shut-down feature known as EcoCalendar. This feature helps minimize utility consumption and HVAC load through a software controlled, automatic steam shut-off program. Utilities can be programmed for automatic start-up and shut-down based on time of day and day of the week.

#### Auto Idle Shut-Off

Like the EcoCalendar, this feature helps save energy by shutting the autoclave off when not in use. It functions similarly to a "sleep-mode" for your home computer. Simply set the desired idle time and the autoclave will enter Sleep Mode after the set idle time has elapsed.

#### EcoJacket™

EcoJacket automatically lowers the temperature of the sterilizer jacket in between cycles, reducing steam and water usage, as well as heat-load to the room.

## Steam Sources

Consolidated sterilizers can operate on house steam or steam produced by integral or remote steam generators. Incoming water quality is always a primary consideration in selection of a steam generator (see Table 9).

Selection of the steam source is based on facility steam availability and other factors associated with sterilizer use and throughput.

### House Steam

In most situations standard house steam, when available, is adequate. An inlet steam filter may be required to clean the steam prior to injection into the sterilizer plumbing system.

### Electric Steam Generator

When house steam is not available, an electric boiler may serve as the steam source. Higher wattage generators produce a higher rate of steam to the chamber. Standard voltages available are 208, 240, 380, 480. Three-phase or single-phase voltage connections are available.

- **Integrated Boiler**  
Consolidated can provide an integral (fits under the sterilizer chamber) steam generator up to 45 kW.  
A standard steam boiler, constructed of carbon steel, is utilized on units fed with tap, or RO water. (See Table 9 for water quality requirements).  
A clean steam boiler, constructed of type 316L stainless steel, is required when the feed water quality is  $>1 \text{ M}\Omega\cdot\text{cm}$ . This option typically requires that the sterilizer be constructed with a stainless-steel chamber, jacket, and plumbing.
- **Remote Boiler**  
Steam generators larger than 45 kW are typically too large to fit beneath the sterilizer chamber and must be remote mounted. Remote boilers are available up to 180 kW.

### Electric/Steam Combination

Sterilizers can be configured to accommodate both house steam and an integral electric generator. This feature allows for easy switching between either steam sources.

### Generator Blowdown

On carbon steel boilers, the generator blowdown function is required to flush impurities from the system and prevent collection of contaminants on the heating elements. The blowdown function can be initiated manually or ordered as an automatic blowdown function if desired.

## Water Quality Compatibility (Stainless Steel vs. Non-Stainless Selection)

Materials that contact the water supply to the sterilizer, including the steam generator, water-cooled devices, water-actuated accessories, valves and piping must be compatible with the purity of the water. See Table 9.

- If deionized or distilled water or water with a purity greater than  $1 \text{ M}\Omega\cdot\text{cm}$  is supplied, then all such points of contact must be of stainless steel construction.
- If water purified by reverse osmosis or tap water with purity of less than  $0.2 \text{ M}\Omega\cdot\text{cm}$  is supplied, then such points of contact must be of carbon steel, bronze or brass construction, as appropriate.

## Air Removal

Depending on the application and required cycle, air removal may be achieved by gravity displacement or mechanical assistance through vacuum for pre- and post-cycle use as required.


- For gravity air removal, the buoyancy of steam is used to displace the air.
- For pre-cycle vacuum, Consolidated offers a choice of a water-ejector vacuum system, or a water-minimizing liquid ring pump.
- For post-cycle vacuum where drying assistance is desired, the sterilizer can use a water-ejector vacuum system or a water-minimizing liquid ring pump.
- Following completion of the post-vacuum function and throughout return to ambient pressure equilibrium, the returning air is filtered through a 0.3 micron bacteria-retentive HEPA filter.

## Shelving

Various shelving options are available upon request. Please see Accessories section on page 9 for further details.

## Compliance and Certifications

Depending on desired application, Consolidated sterilizers can be certified for compliance to the following standards for performance and safety.

- Buy America Act compliant 
- ASME (American Society of Mechanical Engineers): Section VIII, Division I (Unfired Pressure Vessel) code.
- ASME: Section I Code (Miniature Boilers).
- ASME: CSD-1 Code (Controls and Safety for Boilers).
- CRN (Canadian Registration Number): requirements for pressure vessels.
- IEC (International Electrotechnical Commission): EN 61010-1:96—Safety requirements for electrical equipment for measurement, control, and laboratory use. EN 61010-2-041:96—Specific requirements for steam autoclaves.
- UL and cUL (Underwriters Laboratory): Standard 61010-1 and 61010-2.
- US NEC (National Electrical Code) and National Plumbing Code; NFPA-70.
- EN ISO 17665-1: 2024(R2024)—AAMI (Association for the Advancement of Medical Instrumentation) Industry Standard for Moist Heat Sterilization.
- 21 CFR Part 11 (US Code of Federal Regulations, Section 21—Food and Drugs; Electronic Records; Electronic Signatures).
- ANSI/AAMI ST-8: Hospital Steam Sterilizers.

## Consolidated Sterilizer Systems — The X1™ Control System

Consolidated's X1™ controllers offer proven reliability, based on an industrial PLC platform programmed specifically for each Consolidated sterilizer based upon the application. The X1 control system has an intuitive user interface with modern touchscreen display and can be equipped with up to 50 programmable cycles. The entire system is built from industry-standard, open-source components for international availability of parts and service.

Operator Interface	
Display	7" (17.8 mm) Widescreen Color Touchscreen
Resolution	WVGA 800 x 400
Mounting Options	Front or Remote Panel

Industry Standard Sterilizer Cycles	
<ul style="list-style-type: none"> <li>Gravity</li> </ul>	<ul style="list-style-type: none"> <li>Liquid</li> </ul>
Optional Sterilizer Cycles	
<ul style="list-style-type: none"> <li>Pre-Vacuum</li> <li>Liquid with Controlling Load Probe</li> <li>F<sub>0</sub> with Controlling Load Probe</li> <li>Low Temperature (Isothermal)</li> <li>Continuous (Product Lifecycle Testing)</li> <li>Rapid Cool</li> <li>Temperature/Pressure Ramping</li> </ul>	<ul style="list-style-type: none"> <li>Bowie-Dick (Air Removal Test)</li> <li>Vacuum Leak Test</li> <li>Air-Over-Pressure</li> <li>Steam-Air Mix</li> <li>ATF Bioreactor</li> <li>USP 660</li> <li>Customized to Application</li> </ul>

Green Features	
WaterEco® Water Conservation & Monitoring System	Standard
Energy Saver EcoCalendar for Scheduled On/Off Steam Supply	Standard
Auto Idle Shut-Off	Standard
EcoJacket	Standard

Documented Cycle Data Information	
User Defined Cycle Name	Standard
Start Time and Date	Standard
Cycle Type	Standard
Chamber Temperature/Pressures	Standard
Cycle Summary	Standard
Operator ID (If Applicable)	Standard
Cycle Set-Points and Parameters	Standard
Alarms	Standard

Sterilizer Control	
Configurable Cycle Parameters	Standard
Improper Parameter Rejection	Standard
E-Stop	Standard
Three Point Calibration of Sensors	Standard
Automatic Condensate Exhaust Management for all Cycles	Standard
Battery Backup of Program, Memory, & Configuration Data	Standard
Ability to Favorite Cycles	Standard

Security and Process Integrity	
Multi-layered Password Security for Manager and Service	Standard
Ability for Manager to Lock Cycle Configuration	Standard
User Traceability Password Protection	Standard
Up to 50 Usernames and Passwords	Standard

Visual and Audible Alarms	
Cycle Complete, Sensor Failure, Over Temperature, Under Temperature, Over Pressure, Time Limit, Door Seal	Standard
Alarm History Report on Touchscreen	Standard
Preventative Maintenance & Service Reminders	Standard

System Monitoring and Data Storage	
Thermal Printer	Standard
Electronic Data Storage via USB	Standard
Dry Contacts for Door Open, Alarm Condition, and/or Sterilizer On	Optional
SteriNET™ Connex (Remote Troubleshooting) via Ethernet	Optional



## Sterilization Cycles

Consolidated sterilizers can be easily configured for numerous cycles and functions. Each cycle is customizable to meet user specific requirements. All cycle parameters are easily viewed prior to and during the cycle run. The following cycles are pre-configured and installed on all sterilizers:

### Gravity

The gravity cycle uses gravity to displace the air in the chamber with pressurized steam. This cycle is used mainly for unwrapped and nonporous goods. Depending on unit size, the operator may select a sterilization temperature anywhere between 212° F and 275° F\* (100° C and 135° C\*), as well as sterilization time and dry time (see Figure 1).

### Liquid

The liquid cycle prevents boil-over when vented liquids are sterilized. The controller is factory-configured to automatically release steam pressure at an adjustable exhaust rate. Depending on unit size, the operator may select a sterilization temperature anywhere between 212° F and 275° F\* (100° C and 135° C\*), as well as sterilization time (see Figure 2).

Automatic Jacket Blowdown is standard on all sterilizers and is activated during liquid cycles. This feature is used to exhaust steam in the chamber as well as the jacket to prevent liquid loads from boiling over. Factory installed.

### Additional Sterilization Cycles (available upon request)

#### Gravity with Post-Vacuum Drying

This cycle draws a vacuum following sterilization and significantly decreases drying time. Post-vacuum is usually used for wrapped goods, fibers, rubber, bedding, glassware and similar loads. The operator may select a sterilization temperature anywhere between 212° F and 275° F\* (100° C and 135° C\*), as well as sterilization time and dry time (see Figure 3).

#### Liquid with Load Probe

This cycle uses a temperature probe placed within the liquid load to detect, display and print temperature. Applications include sterilization of temperature-sensitive liquids or large volume liquid sterilization (i.e. >2 liter flasks).

#### Pre-Vacuum

This cycle performs a series of steam pulses and vacuum draws prior to sterilization to assure that air has been removed from the load. This also permits the ability to vacuum dry. The number of pulses, pulse steam pressure, vacuum pulse pressure, sterilization time and temperature, and dry time are customizable for each cycle (see Figure 4).

#### Bowie-Dick (Air Removal Test)

Available on pre-vacuum units only. This air-removal test cycle validates the vacuum function using ad hoc test packs as per requirements listed in AAMI ST-8.

#### Vacuum Leak Test

Available on pre-vacuum units only. This cycle provides easy verification of vacuum depth, vacuum seal and chamber piping integrity. Detects leaks greater than 1 mmHg (1 Torr) per minute.

### Effluent Decontamination

Retains liquid effluent, if any, within the chamber for the duration of the sterilization cycle. Aerosol effluent passes through a 0.2 micron hydrophobic HEPA filter with >99.999% efficiency; the filter is located within the chamber and is sterilized in situ to eliminate the need for disposal after each cycle. The filter is easily replaced by the user when required without special tools or service personnel.

### Air-Over-Pressure

This cycle uses compressed air to maintain chamber pressure at or above the sterilization pressure until the liquid has cooled to a user adjustable temperature. This is useful for small volumes of liquid susceptible to boil-over (see Figure 5). Air is provided using a house air source.

### Low Temperature

Typically used for temperature-sensitive goods, this function creates a zero-pressure, temperature only cycle between 190° F and 212° F (88° C and 100° C). Often referred to as an isothermal or inspissation cycle.

### Rapid Cooling

Useful for rapidly cooling liquid media via an interior water spray mechanism or jacket quench, post-sterilization.

### F<sub>0</sub> Cycle with Load Probe

Useful for heat-sensitive liquid media. Sterilization begins when temperature reaches 212°F and is completed when the desired F<sub>0</sub> set-point is reached. F<sub>0</sub> is adjustable.

### Temperature/Pressure Ramping

Allows the user to specify a) the amount of time it takes for the load to rise to sterilization temperature/pressure and b) the amount of time it takes to exhaust the sterilizer chamber post-cycle. Ideal for loads that are sensitive to rapid pressure changes.

### Continuous/Product Life Cycle Testing

Any sterilization cycle can be programmed to be continuously repeated (number of times is adjustable) without any input from the operator. Used for product lifecycle testing.

### Steam-Air Mix

Useful for liquid-filled syringes or other sealed containers. Sterilization occurs at an elevated pressure relative to temperature. Extra pressure is achieved using a house air source.

### ATF Bioreactor Cycle

The ATF bioreactor cycle is designed to permit the sterilization of hollow fiber filters while housed in bioreactors. ATF (Alternating Tangential Flow) is a technology used in the perfusion culture of mammalian cells, proteins and antibodies. This cycle incorporates a series of temperature ramps and hold times to prevent damage to the fiber filter.

### USP 660 Cycle

The USP 660 cycle is used for sterilizing glass containers for pharmaceutical packaging as defined in the US Pharmacopeia in Chapter <600> Containers. Comes with load probe.

### Custom Cycles

Custom cycles are available upon request. Please contact Consolidated to further discuss your application.

\* Exact upper temperature limit is model specific and may vary depending on model and/or options. Contact Consolidated for a solution tailored to your requirements.

## Sterilizer Cycle Selection

### Gravity. (Figure 1)

Steam flows through the sterilizer; temperature and pressure ramps slightly to a continuous flow purge.

The controller closes the exhaust valve and interior temperature and pressure ramp up to desired setpoint.

The program maintains dwell until desired time is reached, when pressure is released.

The interior is restored to ambient pressure although contents remain relatively hot.

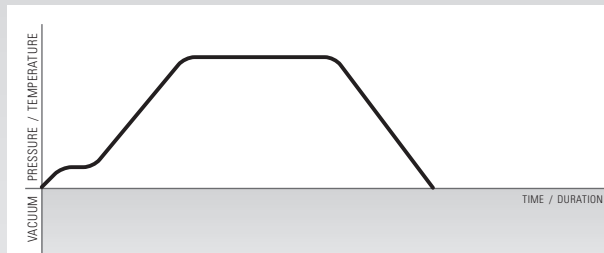


Figure 1: Gravity Cycle

### Liquid. (Figure 2)

Steam flows through the sterilizer; temperature and pressure ramps slightly to a continuous flow purge.

The controller closes the exhaust valve and interior temperature and pressure ramp up to desired setpoint.

The program maintains dwell until desired time is reached. Pressure is gradually released to allow dissolved gases to reach equilibrium to prevent boil-over.

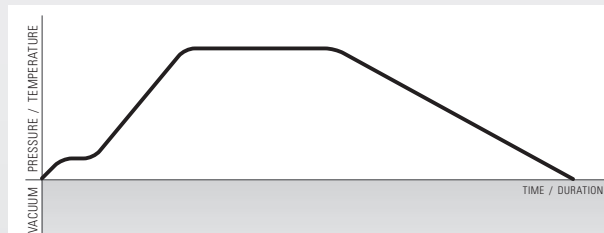


Figure 2: Liquid Cycle

### Gravity with Post-Vacuum Drying. (Figure 3)

Steam flows through the sterilizer; temperature and pressure ramps slightly to a continuous flow purge.

The controller closes the exhaust valve and interior temperature and pressure ramp up to desired setpoint.

The program maintains dwell until desired time is reached, when pressure is released.

A vacuum system draws the interior atmosphere through ambient to negative pressure, accelerating drying time as programmed.

Once completed the vacuum is released and the load is restored to ambient pressure.

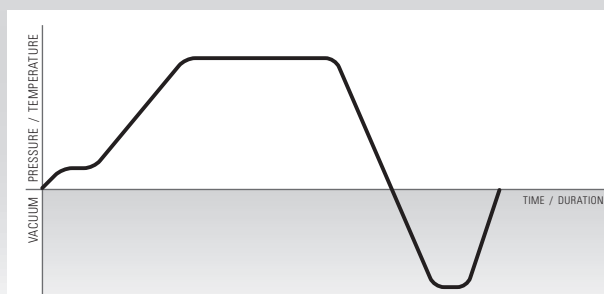


Figure 3: Gravity with Post-Vacuum Drying Cycle

### Pre-Vacuum with Post-Vacuum Drying. (Figure 4)

Steam flows through the sterilizer; temperature and pressure ramps slightly to a continuous flow purge.

The controller closes the exhaust valve and interior temperature and pressure ramp to an intermediate setpoint without dwell.

A vacuum system draws the interior atmosphere to negative pressure to remove latent air from concealed pockets within the load.

The process is repeated per program and protocol. Following the final pre-vacuum pulse, the sterilizer ramps to desired setpoint.

The program maintains dwell until desired time is reached, when pressure is released.

If a post-vacuum program is used, the vacuum system draws the interior atmosphere through ambient to negative pressure, accelerating drying time as programmed, after which the vacuum is released and the load is restored to ambient pressure.

If a post-vacuum program is not used, the load is restored to ambient pressure.

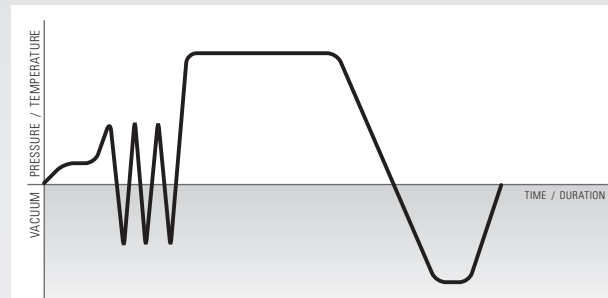


Figure 4: Pre-Vacuum with Post-Vacuum Drying Cycle

### Liquid, Air-Over-Pressure. (Figure 5)

Usually used with smaller amounts of media to prevent liquid loss.

The air-over-pressure function operates identical to the liquid cycle except that pressurized air is injected to cool the load while maintaining pressure to displace steam.

After sterilization, air is injected, pressure is maintained, this prevents boil-over and minimizes evaporation.

Air pressure is then released and the load is restored to ambient pressure.

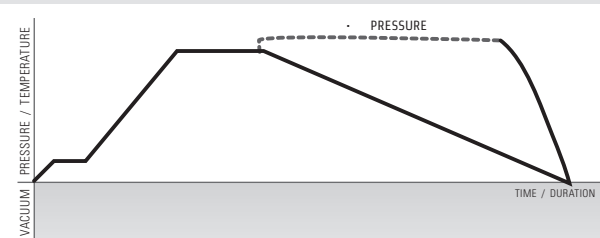


Figure 5: Liquid, Air-Over-Pressure Cycle

## Options

Consolidated offers a range of options and accessories to optimize and customize your sterilizer, to meet requirements for internal or third-party compliance, and to improve workflow. Options are factory installed and must be specified when ordering. For further information about available options and tailoring your sterilizer to meet your needs contact Consolidated.

### Sterilizer Configuration Options

#### Chamber Finish

Chambers can be provided with a passivated and electropolished and/or high polished to less than 10Ra interior surface.

#### Nickel-Clad Chamber

Recommended chamber material for sterilizing high-concentration salt solutions or if the feed-water or tap-water for steam generation has high chloride content. Nickel-Clad chambers are provided with a carbon steel jacket.

#### Interior Chamber Tracks

Chamber tracks may be installed into the sterilizer to permit use of an existing CSS cart & carriage or use with a future cart and carriage. Chamber tracks must be factory installed.

#### Controller Mounting

Choose where the sterilizer control system is mounted: left-side, right-side or remote-mounted on an adjacent wall.

#### Plumbing Side

Depending on the area surrounding the autoclave, and to make it easy for your service team, choose the desired location (left or right side) of the sterilizer's components and piping.

#### Stainless Steel Piping

Upgrade from brass/bronze piping & components to type 316 stainless steel for all product wetted surfaces.

#### Other Options:


Add your organization's logo to the autoclave, add additional shelving, add extra sampling ports, request a custom frame width, and many more!

### Generator Options

#### Automatic Generator Blowdown

Recommended to remove accumulated impurities from the carbon steel with a steam generator.

### Water and Energy Saving Options

 **Steam Sterilizer Water Conservation & Monitoring Systems** The Consolidated WaterEco® Systems reduce water consumption to the autoclave by up to 99%. All WaterEco® systems communicate with the sterilizer control system to provide real-time data for cooling water consumed per cycle, day, month, and year. The system also triggers an alarm (displayed on the touchscreen, printout, and history report) if the waste temperature exceeds set-point or if there is a sensor failure. Factory or field installed. Available on gravity and pre-vacuum units in the following configurations:

- **WaterEco®:** This device is included on all Consolidated sterilizers and reduces cooling water usage by up to 90%. In addition to saving water, it monitors water consumption through the touchscreen display. The system is designed to cool effluent using a minimal amount of cold water, air, and previously cooled effluent.
- **WaterEco® Plus:** Reduces cooling water consumption by up to 99% through a stainless steel heat exchanger that utilizes a facility chilled water supply. This system virtually eliminates once-through cooling water.
- **WaterEco® Vac Plus:** This full recovery system reduces water consumption by the vacuum and cooling systems by up to 99%. Designed to integrate with pre-vacuum autoclaves when facility chilled water is readily available. This system is the best solution to minimize water consumption.

#### Jacket Steam Return

Sterilizer is configured with a jacket steam return connection port. This helps save energy and reduce steam consumption by returning sterile steam from the jacket back to the "house" facility boiler through a user-supplied condensate return line. For house steam units only.

#### Form C Dry Contacts — Hood Control

This feature integrates the sterilizer with a local canopy hood to reduce unnecessary venting of conditioned air. Form C Dry Contacts signal your building automation system to power-down the hood when the door of the autoclave is closed.

### System and Data Monitoring



#### **SteriNET® Remote Monitoring and Troubleshooting**

SteriNET® Connex permits Consolidated engineers to provide quick, accurate and economical offsite diagnosis of operation problems through an Ethernet connection. This option is ideal for facilities located in remote areas where service assistance is hours or days away. This option includes one year of free, remote-transmitted software upgrades from the Consolidated technical support center.

#### **Form C Dry Contacts**

Offers dry connection to building monitoring system. Specify desired outputs when ordering: door open/close, power on/off or sterilize alarm conditions.

### Probe and Access Port Options

#### **Load Probe**

Includes a temperature probe with lead wire permitting placement in the load to be sterilized such as liquid, red-bag or other. Sterilization cycle time is based on load temperature instead of chamber temperature. Factory installed.

#### **Multi-Probe Sealing Gland**

Multi-port gland permits safe installation of thermocouples or other probes inside the chamber. Factory installed. Specify desired quantity and location when ordering.

#### **Validation Port**

All CSS sterilizers are equipped standard with one 1/2" NPT opening into the chamber. An additional port of up to 1" NPT or 2" tri-clamp may be added.

### Safety Options

#### **Flood Switch**

In the event of excessive liquid in the chamber or jacket, an alarm will sound and a message will be displayed warning the user about the flood condition. Factory installed.



### Seismic Anchorage Details

May be required to meet building codes in seismically active areas. Contact your facility manager to determine if required. If ordered, Consolidated will provide anchor attachment points and suggested seismic anchorage details.

### Pass-Thru (Double-Door) Options

#### Dual Controls

For pass-thru models only. Provides full touchscreen controls mounted on each end of the unit for operation from within or outside the containment area.

#### Pass-thru Door Interlock System

For pass-thru models only. Prevents opening of the sterilizer door on the non-contained side until completion of a sterilization cycle. Also prevents both doors from being opened simultaneously. This system is programmed to accommodate the desired direction of product flow and is equipped with a secondary touchscreen controller located on the non-contained side.

#### Effluent Decontamination Control

Exclusively from Consolidated, this effluent decontamination system traps bacteria within the autoclave chamber and sterilizes it prior to release into the drain (Consolidated Patented Process). Any liquid effluent is held in the chamber for the duration of the sterilization cycle. All aerosol effluent passes through a 0.2 micron hydrophobic HEPA filter with >99.999% efficiency. The steam that sterilizes the products in the chamber also sterilizes the bacteria held by the filter. Filters are easily changed after use. Factory installed. Required for BSL-3 applications.

#### Bioseal® Biological Sealing Flange

The Bioseal® biological sealing flange, required for Biological Safety Level-3 (BSL-3) applications, employs a stainless steel barrier flange welded to the periphery of the sterilizer and another stainless steel barrier attached to the wall opening through which the sterilizer is installed. This arrangement is used to prevent the passage of airborne microorganisms from the contained side of the sterilizer to the uncontained side. A single piece solid-silicone gasket completes the seal between the building and the sterilizer thereby ensuring seal integrity is not compromised during normal expansion and contraction of the building.

#### Air Differential Seal

An Air Differential Seal is required for applications that require a negative or positive air pressure differential between two rooms adjoined with an autoclave. The seal prevents the bulk passage of air between areas maintained at a pressure differential. A stainless steel barrier flange is welded to the periphery of the sterilizer and solid-silicone gasket secures the welded seal to the wall and permits normal expansion and contraction of the building.

#### Vermin Seal

A vermin seal employs a metal barrier to impede rodents from passing through the space between a recessed-mounted autoclave and the wall opening. This is useful if a recessed autoclave is located in the same room as live rodents.

#### BSL-3 Package

Required options for pass-thru autoclaves installed in Biosafety Level 3 (biocontainment) labs:

- Dual Controls
- Pass-thru Door Interlock System
- Effluent Decontamination
- Bioseal
- Up to 50 Password Protected Usernames

- Lockable Cycle Parameters

For more information on BSL-3 configurations, options, and requirements see the Guide to BSL3 Autoclaves E-Book.

### Accessories

#### Water Purification System

High-purity water purification systems are available for facilities that do not meet the minimum water purity requirements (see Table 9) or that require clean steam. Specify when ordering.

- **Option 1:** A reverse osmosis (RO) water filtration system designed to pre-treat autoclave feed-water can be supplied (Model WPS-1200-RO).
- **Option 2:** If deionized water is required for clean steam generation, an RO/DI system that incorporates extra filtration to produce Type II deionized water (>1 MΩ resistivity) can be supplied (Model WPS-1200-DI).

#### Shelving

Shelves and racks are constructed of electro-polished type 316L stainless steel. Specify when ordering.

- **Option 1:** One (1) stationary full-length wire-grid bottom shelf and one (1) upper stationary shelf suspended in the middle of the chamber on shelf brackets.
- **Option 2:** One (1) stationary full-length wire-grid bottom shelf and two (2) upper stationary shelves suspended on two evenly spaced shelf brackets.
- **Option 3:** One (1) extendable full-length wire-grid bottom shelf and one (1) upper extendable shelf suspended in the middle of the chamber on shelf brackets.
- **Option 4:** One (1) extendable full-length wire-grid bottom shelf and two (2) upper extendable shelves suspended on two evenly spaced shelf brackets.

#### Loading Cart & Transfer Carriage

A loading cart, mounted on rollers, with two stainless steel shelves and a wheeled transfer carriage. The transfer carriage mates to guides on the interior chamber tracks to ease loading and unloading of sterilization loads up to 500 lbs evenly distributed. Chamber tracks must be configured in chamber at time of build to utilize loading cart and carriage; specify when ordering.

#### Uninterruptible Power Supply (UPS)

Allows the control system to remain on in the event of a brief power loss or brownout. Specify when ordering.

#### Sump Pump

Required when gravity drain is not available. Provides positive pump to remove condensate effluent to hard-plumbed drain. Specify when ordering.

#### Steam Inlet Filter

Recommended to clean dirty steam (i.e. remove particulates) from facility steam sources. Specify when ordering.



## Validation

Validation protocols and/or execution of protocols is available from Consolidated for new or existing autoclaves. Steril-Q™ Validation Services & Documentation includes an array of documentation, tools and expertise required to successfully qualify and validate an autoclave within accepted protocols that comply with the current USA and international standards for steam sterilization as set forth in ISO 17665. The various protocols and services include:

- **Installation Qualification (IQ):** Provides documented evidence that the equipment has been built and installed to specification, and that all supporting services (i.e., utilities such as electricity, water, and steam) are available and connected properly.
- **Operation Qualification (OQ):** Provides documented evidence that the sterilizer operates in accordance with design specifications.
- **Performance Qualification (PQ):** Documents the sterilizer's ability to achieve the desired outcome (i.e. sterilization of specific loads when operated in accordance with pre-defined operating procedures).
- **Steam Quality Testing (SQ):** Documents the suitability of the steam delivered to the chamber, ensuring compliance with the requirements of ISO 17665. A successful test result ensures a predictable level of process lethality.
- **Factory Acceptance Testing (FAT):** Evaluates the sterilizer after assembly to verify that its operation complies with design specifications prior to shipment from Consolidated's corporate headquarters.
- Onsite validation and execution of protocols as well as cycle development services are available. For more information see the Autoclave Validation E-Book.

## Site Preparation and Utilities

### General

Utilities of ample capacity required for operation of the sterilizer shall be provided at the place of installation at the facility. Key considerations include electric power availability, steam source, water type and flow capacity, and compressed air (if air-over-pressure or steam-air mix cycle is desired). For information on required utilities and locations refer to the sterilizer architectural drawings and Tables 3–6, 9.

- Drain and ventilation considerations include wastewater and the facility HVAC system.
- Clearances include door and cart allowances, recommended 18–24" for access to plumbing, adequate load-bearing capacity of floor at installation site, and biointegrity.
- Consider the location of sprinklers, smoke detectors, and canopy capture hoods.
- Contact Consolidated or your Consolidated sales representative for planning and selection assistance.

## Installation

Consolidated offers a choice of installation options to accommodate the facility schedule. Permitting and inspections are not included. In general, Consolidated can arrange for any of the following installation preferences:

- **Level 1:** Sterilizer start-up and user training.
- **Level 2:** Sterilizer uncrating, setting in place, leveling, final assembly, start-up, and user training.
- **Level 3A:** Sterilizer uncrating, setting in place, leveling, final assembly, final utility connections, start-up, and user training.
- **Level 3B:** Receiving the shipment, sterilizer uncrating, setting in place, leveling, final assembly, start-up, and user training.
- **Level 4 (Turnkey):** Receiving the sterilizer, uncrating, setting in place, leveling, final assembly, final utility connections, start-up, and user training.
- **Installation Supervision:** Consolidated will supervise on-site installation work performed by facility designated labor.

## Preventative Maintenance

At Consolidated, we believe that sterilizers should be easy to use, service, and maintain. To maximize uptime and keep your equipment in peak condition, we offer SterilCARE™ service plans and PartsPAX solutions to simplify autoclave maintenance.

### Parts


Consolidated offers the convenience of three PartsPAX for both the steam sterilizer and integral steam generator. Each PartsPAX includes the recommended certified components to keep the sterilizer working reliably and safely over time.

- **Safety PAX:** Includes components that ensure your autoclave is running safely. Recommended to be changed out annually regardless of number of cycles.
- **Performance PAX:** Includes all essential components to ensure the sterilizer is running smoothly. Recommended to be changed out every 300 cycles.
- **Complete PAX:** Includes all Performance PAX and Safety PAX components plus a few other key recommended replacement parts. Recommended to be used if maintenance hasn't been performed in over a year and/or the autoclaves sees heavy use.

### Service

Consolidated offers two simple SterilCare service plans: Safety Plus and Proactive. Consolidated Sterilizer Systems recommends a safety inspection annually (Safety Plus) and a preventative maintenance inspection by an authorized factory trained service representative once every 300 cycles (Proactive).

### Standard Mounting Configurations - Left Hinge, Right Controls

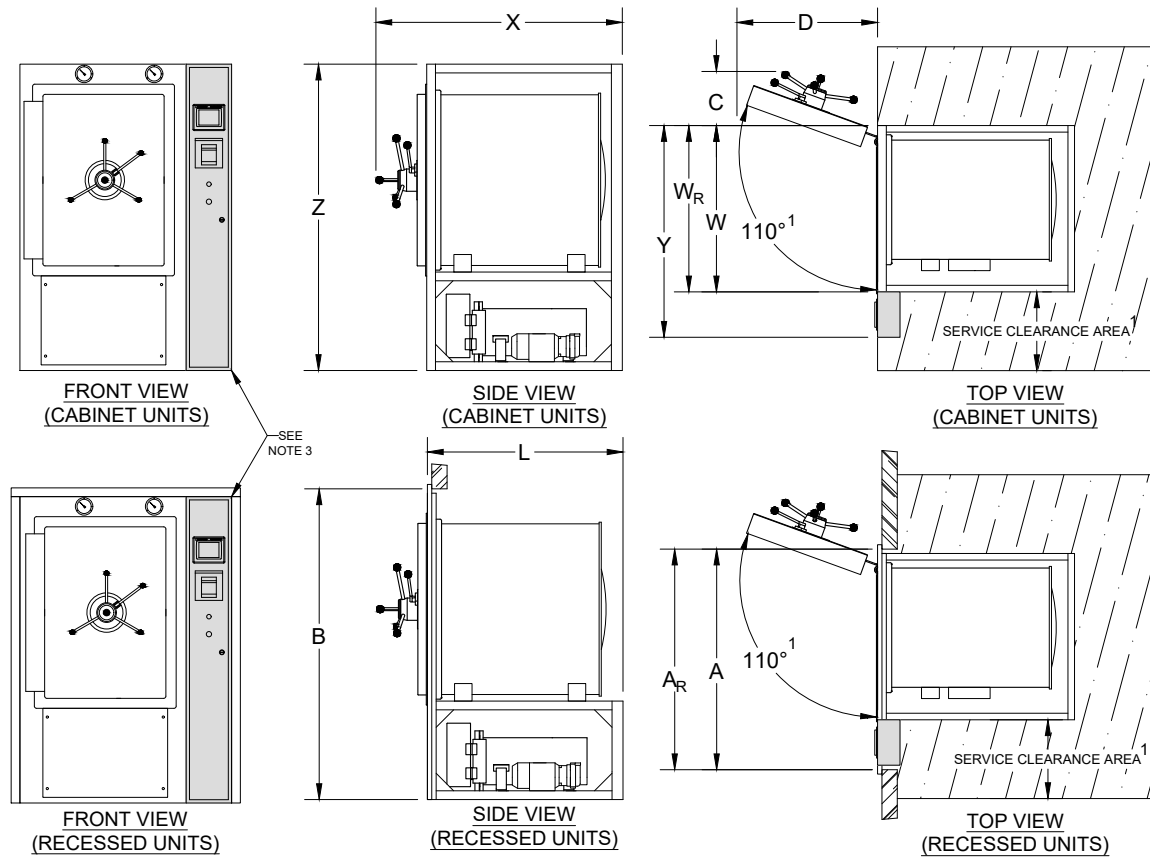

**Table 1: Sterilizer Dimensions<sup>2</sup>**

Model	SR-24C	SR-24D	SR-24E <sup>6</sup>	SR-24F <sup>6</sup>	SR-26B
Chamber Dimensions (w x h x f-b)	24" x 36" x 36" 61 x 91.4 x 91.4 cm	24" x 36" x 48" 61 x 91.4 x 122 cm	24" x 36" x 60" 61 x 91.4 x 152.4 cm	24" x 36" x 72" 61 x 91.4 x 183 cm	26" x 26" x 49" 66 x 66 x 124.5 cm
Volume	18 cu. ft (510L)	24 cu. ft (680L)	30 cu. ft (850L)	36 cu. ft (1019L)	19.2 cu. ft (544L)
Overall Length (X)	56" 142.2 cm	68" 172.7 cm	80" 203.2 cm	92" 233.68 cm	68" 172.7 cm
Overall Width (Y)	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm
Overall Height (Z) <sup>5</sup>	71" 180.3 cm	70.25" 178.4 cm	70.25" 178.4 cm	70.25" 178.4 cm	71" 180.3 cm
Frame Length (L)	45" 114.3 cm	58" 147.3 cm	70" 177.8 cm	82" 208.3 cm	55" 114.3 cm
Frame Width (W)	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm
Frame Width - Restricted (W <sub>R</sub> ) <sup>7</sup>	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm
Wall Opening Width (A)	50.375" 128 cm	50.375" 128 cm	50.375" 128 cm	50.375" 128 cm	50.375" 128 cm
Wall Opening Width - Restricted (A <sub>R</sub> ) <sup>7</sup>	53.375" 135.6 cm	53.375" 135.6 cm	53.375" 135.6 cm	53.375" 135.6 cm	53.375" 135.6 cm
Wall Opening Height (B)	72" 182.9 cm	71.25" 181 cm	71.25" 181 cm	71.25" 181 cm	72" 182.9 cm
Door Swing Clearance (C)	20" 50.8 cm	20" 50.8 cm	20" 50.8 cm	20" 50.8 cm	12.75" 32.4 cm
Door Swing (D)	31.5" 80 cm	31.5" 80 cm	31.5" 80 cm	31.5" 80 cm	35.5" 90.2 cm

1. Recommended service clearance is 18-24" both sides and back. If necessary, service clearance can be decreased or adjusted to one side to accommodate facility space constraints.
2. Additional options may require a larger footprint.
3. The control housing is shipped detached from the sterilizer to allow passage through doorways, reducing pre-installation Overall Width (Y) by 10.375". When the sterilizer is installed, the control housing and electrical connections are easily attached.
4. All views contain configuration specific components. These are for illustrative purposes only, actual configuration may vary.
5. After adjustment, leveling feet may add up to 1 inch to the Overall Height (Z).
6. Electric generator is not integral to sterilizer and requires a separate footprint.
7. Restricted frame and wall opening is required when there is no service on the hinge side. Contact CSS for more info.
8. Left hinge and right controls is shown. Opposite hinge and controller side is available.

This page shows a variation of recessed mounting to allow a sterilizer to fit into an existing narrow wall opening

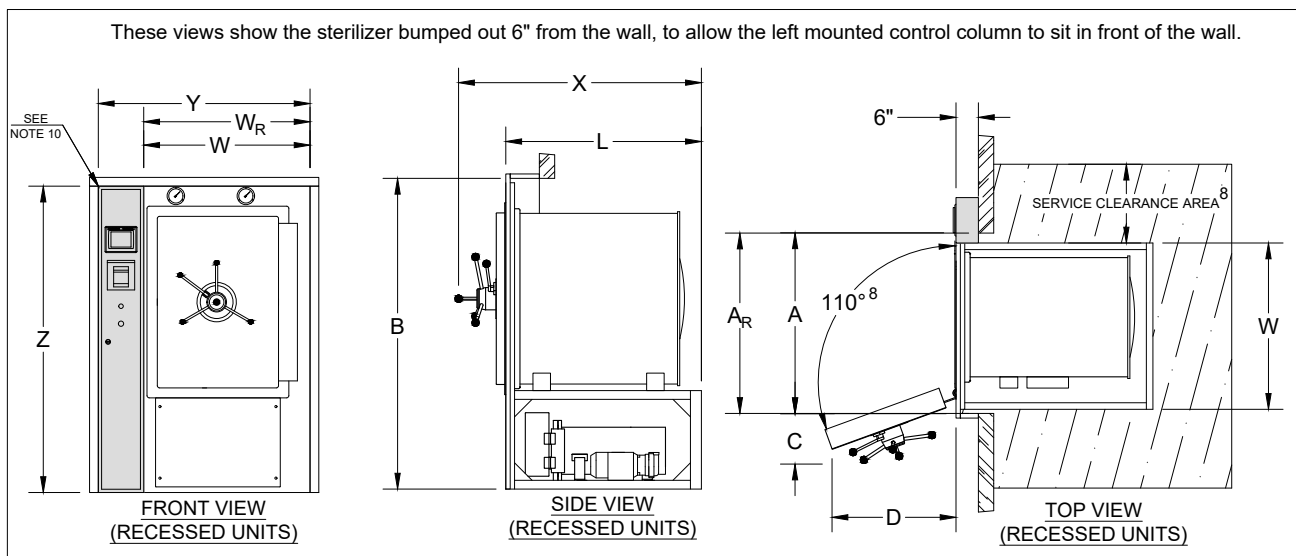


Table 2: Sterilizer Dimensions<sup>9</sup>

Model	SR-24C	SR-24D	SR-24E <sup>11</sup>	SR-24F <sup>11</sup>	SR-26B
Chamber Dimensions (w x h x f-b)	24" x 36" x 36" 61 x 91.4 x 91.4 cm	24" x 36" x 48" 61 x 91.4 x 122 cm	24" x 36" x 60" 61 x 91.4 x 152.4 cm	24" x 36" x 72" 61 x 91.4 x 183 cm	26" x 26" x 49" 66 x 66 x 124.5 cm
Volume	18 cu. ft (510L)	24 cu. ft (680L)	30 cu. ft (850L)	36 cu. ft (1019L)	19.2 cu. ft (544L)
Overall Length (X)	56" 142.2 cm	68" 172.7 cm	80" 203.2 cm	92" 233.68 cm	68" 172.7 cm
Overall Width (Y)	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm	48.375" 122.9 cm
Overall Height (Z)	71" 180.3 cm	70.25" 178.4 cm	70.25" 178.4 cm	70.25" 178.4 cm	71" 180.3 cm
Frame Length (L)	45" 114.3 cm	58" 147.3 cm	70" 177.8 cm	82" 208.3 cm	55" 114.3 cm
Frame Width (W)	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm	38" 96.5 cm
Frame Width - Restricted (W <sub>R</sub> ) <sup>12</sup>	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm	41" 104.1 cm
Wall Opening Width (A)	39" 99.2 cm	39" 99.2 cm	39" 99.2 cm	39" 99.2 cm	39" 99.2 cm
Wall Opening Width - Restricted (A <sub>R</sub> ) <sup>12</sup>	42" 106.8 cm	42" 106.8 cm	42" 106.8 cm	42" 106.8 cm	42" 106.8 cm
Wall Opening Height (B)	72" 182.9 cm	71.25" 181 cm	71.25" 181 cm	71.25" 181 cm	72" 182.9 cm
Door Swing Clearance (C)	20" 50.8 cm	20" 50.8 cm	20" 50.8 cm	20" 50.8 cm	12.75" 32.4 cm
Door Swing (D)	31.5" 80 cm	31.5" 80 cm	31.5" 80 cm	31.5" 80 cm	35.5" 90.2 cm

8. Recommended service clearance is 18-24" both sides and back. If necessary, service clearance can be decreased or adjusted to one side to accommodate facility specific space constraints.


9. Additional options may require a larger footprint.

10. The control housing is shipped detached from the sterilizer to allow passage through doorways, reducing pre-installation Overall Width (Y) by 10.375". When the sterilizer is installed, the control housing and electrical connections are easily attached.


11. Electric generator is not integral to sterilizer and requires a separate footprint.

12. Restricted frame and wall opening is required when there is no service on the hinge side. Contact CSS for more info.

## Typical Utility Requirements

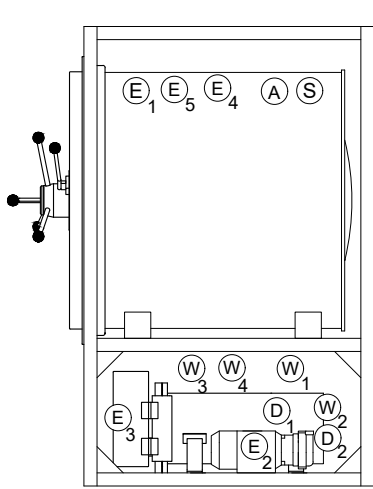

**SIDE VIEW**

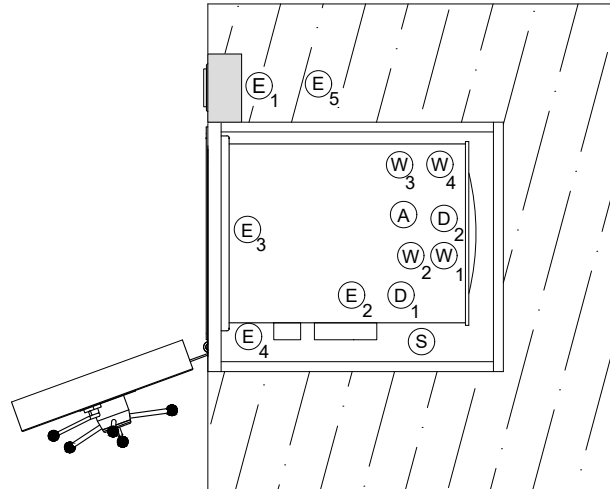

**TOP VIEW**

### UTILITY LEGEND

- (E)** Electrical
- (D)** Drain
- (W)** Water
- (S)** Steam
- (A)** Air

**Table 3: General Connections**

Name	Symbol	Connection on Sterilizer	MEP Requirements <sup>15</sup>
Water Eco/Vacuum System	(W) <sub>1</sub>	1/2" NPT	45 psig dynamic min, 80 psig static max, 3/4"-1" NPT, 12 gpm capacity, with shut off valve and union <sup>13</sup>
Controls System	(E) <sub>1</sub>	NEMA 15-5 Plug	120VAC 10A non-GFCI outlet/220VAC 10A hard wire
Sterilizer Drain	(D) <sub>1</sub>	1 1/2" NPT	2 1/2" air gap, 3" diameter minimum funnel required, 15 gpm, min 1 1/2" drain piping, 6" floor sink recommended, location external to unit footprint recommended
Ethernet Port (Optional)	(E) <sub>5</sub>	RJ-45	Standard RJ-45 ethernet cable with internet access
Direct Steam (Optional)	(S)	3/4" NPT	50-80 psig dynamic, 3/4"-1" NPT, 180 lbs/hr capacity, insulated line with drip leg, shut-off valve and union
Dry Contacts (Optional)	(E) <sub>4</sub>	2-Wire Terminal	2 signal wires per contact compatible with Form C Dry Contacts
Compressed Air (Optional)	(A)	1/2" NPT	50-75 psig, 10-35 SCFM, 99% dry & oil free, backflow preventer not provided by CSS

14. (W)<sub>1</sub> cold water connection will be a single source to feed both water eco system and any vacuum system ordered with the unit. Water quality must comply with the General Vacuum Device & Quench specification in Table 9. Backflow preventer required, not provided by CSS.

**Table 4: Optional Vacuum Systems (maximum one per unit)**

Name	Symbol	Connection on Sterilizer	MEP Requirements <sup>15</sup>
Booster Pump	(E) <sub>2</sub>	Hard Wire	120VAC/208-230VAC, single phase, minimum 20 Amp circuit required
Vacuum Pump - Electrical	(E) <sub>2</sub>	Hard Wire	208/240/480VAC, 3-phase, minimum 20 Amp circuit required

**Table 5: Optional Electric Steam Generator**

Name	Symbol	Connection on Sterilizer	MEP Requirements <sup>15</sup>
Power Supply	(E) <sub>3</sub>	Hard Wire	Available in 208/240/380/480VAC, 3-phase, see Table 7 for amp draw, fused disconnect required, see local codes for remote shut down switch requirements.
Generator Water Feed	(W) <sub>2</sub>	1/2" NPT	45 psig dynamic min, 80 psig static max, 1/2" NPT, 1 gpm, with shut off valve and union <sup>14</sup>
Manual Generator Drain	(D) <sub>2</sub>	1/2" NPT	See D1 in Table 3

15. Generator feedwater must comply with the Generator Water Feed specification in Table 9. Backflow preventer not provided by CSS.

**Table 6: Optional Chilled Water (WaterEco Plus and Vac Plus Units Only)**

Name	Symbol	Connection on Sterilizer	MEP Requirements <sup>15</sup>
Chilled Water Feed	(W) <sub>3</sub>	1" NPT	20 psig dynamic min, 35°F-55°F, 10 gpm insulated with ball valve and union
Chilled Water Return	(W) <sub>4</sub>	1" NPT	-5 psi drop on return side, +15°F max rise

16. For a more detailed list of MEP requirements and specifications see document 92006-01 "Autoclave Installation and Utility Overview".



Table 7: Power and Steam Usage<sup>16</sup>

Power and Steam Usage			Electrically Heated				Steam Heated			
Model	Chamber Dimensions (w x h x f-b)	Air Removal Method	Generator Size (KW)	Generator Current (amps) <sup>17</sup>				Steam Consumption		
				208V	240V	380V	480V	Peak (lb/hr)	Per Cycle (lb/cycle)	Idle (lb/hr)
SR-24C	24" X 36" X 36"	Gravity	30	83	72	46	36	180	40	9
	61 X 91.4 X 91.4 cm	Vacuum	45	125	108	68	54	180	65	9
SR-24D	24" x 36" x 48"	Gravity	30	83	72	46	36	180	45	10
	61 x 91.4 x 122 cm	Vacuum	45	125	108	68	54	180	75	10
SR-24E	24" x 36" x 60"	Gravity	60	167	144	91	72	180	50	11
	61 x 91.4 x 152.4 cm	Vacuum	80-100	222-278	192-240	122-152	96-120	180	105	11
SR-24F	24" x 36" x 72"	Gravity	80	222	198	122	96	180	55	11
	61 x 91.4 x 183 cm	Vacuum	100-120	278-333	240-289	152-182	120-144	180	125	11
SR-26B	26" x 26" x 49"	Gravity	30	83	72	46	36	180	40	9
	66 x 66 x 124.5 cm	Vacuum	45	125	108	68	54	180	70	9

17. Assuming 30 Minute sterilizing time at 250°F (121°C) and 5 minute drying time.

18. Nominal current drawn by a 3-phase generator. Local codes and regulations may affect breaker size. Single phase available if required.

Table 8: Water Consumption (Per Chamber)<sup>18</sup>

Model	Chamber Dimensions (w x h x f-b)	Air Removal Method	Water Consumption					
			Cold Water			Hot/Treated Water		
			Peak (gpm)	Per Cycle (gal/cycle)	Idle (gph)	Peak (gpm)	Per Cycle (gal/cycle)	Idle (gph)
SR-24C	24" X 36" X 36" 61 X 91.4 X 91.4 cm	Gravity	6	5	1	1	5	1
		Ejector	6	80	1	1	5	1
		Vac Pump	6	37	1	1	5	1
SR-24D	24" x 36" x 48" 61 x 91.4 x 122 cm	Gravity	6	6	1	1	5	1
		Ejector	6	97	1	1	5	1
		Vac Pump	6	47	1	1	5	1
SR-24E	24" x 36" x 60" 61 x 91.4 x 152.4 cm	Gravity	6	8	1	5	6	1
		Ejector	6	113	1	5	6	1
		Vac Pump	6	53	1	5	6	1
SR-24F	24" x 36" x 72" 61 x 91.4 x 183 cm	Gravity	6	9	1	5	7	1
		Ejector	6	130	1	5	7	1
		Vac Pump	6	67	1	5	7	1
SR-26B	26" x 26" x 49" 66 x 66 x 124.5 cm	Gravity	6	5	1	1	5	1
		Ejector	6	84	1	1	6	1
		Vac Pump	6	39	1	1	6	1

19. Assuming 30 minute sterilizing time at 250°F (121°C) and 5 minute drying time.

Table 9: Nominal Water Quality Requirements

1 mg/L = 1 ppm	Steam Generator Boiler Feed Water <sup>19, 20, 21</sup>				General Vacuum Device & Quench	
	Boiler Feed Water		Feed Water into CSS RO Filter System		--	--
Characteristic	Recommended Condition	Maximum Condition	Recommended Condition	Maximum Condition	Recommended Condition	Maximum Condition
Temperature [°F(°C)]	As Supplied	140 (60)	-	100 (37.8)	40-60 (4-16)	70 (21)
Silica (ppm)	0.0-1.0	2.5	2.5-15	15	0.0-1.0	2.5
TDS (ppm)	50-150	250	100-1500	1500	50-200	500
pH	7.5-8.5	7.5-9.0	7.5-8.5	7.5-9.0	6.8-7.5	6.5-9.0
Alkalinity (ppm)	50-180	350	50-180	350	50-180	350
Hardness (ppm)	5-25	85	25-225	225	10-85	171
Resistivity (Ω-cm)	2,000	26,000 <sup>27</sup>	-	-	2,000-500,000	500,000
Specific Conductance (µmΩ/cm)	500	38.5	-	-	500-2	2
Sodium (ppm)	0.0-2.5	10	2.5-100	100	-	-
Chloride (ppm)	0.0-2.5	10	2.5-125	125	-	-
Copper (ppm)	0.0-0.2	-	-	-	-	-
Iron (ppm)	0.0-0.2	-	-	-	-	-
Zinc (ppm)	0.0-0.2	-	-	-	-	-
Aluminum (ppm)	0.0-0.2	-	-	-	-	-
Magnesium (ppm)	0.0-0.2	-	-	-	-	-

20. Stainless-steel generators require deionized water with resistivity ≥ 1MΩ•cm.

21. If water supplied is greater than 26,000 Ω•cm contact Consolidated for recommendation.

22. Boiler feed water quality applies to both Integrated and House Steam boilers.

Table 10: Weight & Heat Loss Data			Peak Heat Loss (BTU/hr at 70°F [21°C])		
			Cabinet	Recessed	
Model	Heat Source Options	Max. Operating Weight <sup>21</sup>	To Room	Front of Wall	Back of Wall
SR-24C	Steam	3000 lbs	12550	5000	7500
	Electric	3175 lbs	14800	5800	8950
SR-24D	Steam	3525 lbs	14650	5000	9600
	Electric	3700 lbs	16900	5800	11050
SR-24E	Steam	4075 lbs	16800	5000	11800
	Electric	4075 lbs <sup>22</sup>	19050	5800	13250
SR-24F	Steam	4875 lbs	18950	5000	14000
	Electric	4875 lbs <sup>22</sup>	21200	5800	15450
SR-26B	Steam	2950 lbs	14390	5000	9400
	Electric	3125 lbs	16140	5700	10500

23. Assuming chamber fully loaded with flasks filled 25% with water.

24. Remote Mounted Generator.

[illegible]